



# Instructions for Use

380349-B

English

For Research Use Only

## RUO LZ-SAA Calibrator EIKEN

**REF** V-SZ12

### INTENDED USE

RUO LZ-SAA Calibrator EIKEN is intended for use with RUO LZ TEST EIKEN SAA (**REF** V-SZ11), RUO QC-SAA L EIKEN (**REF** V-SZ13) and RUO QC-SAA H EIKEN (**REF** V-SZ14) as the calibrator when using automated analyzers to measure Serum Amyloid A (SAA) in serum or plasma. For research use only, not for use in diagnostic procedures.

### CALIBRATOR VALUE

The calibrator material is standardized against the WHO International Standard (NIBSC code : 92/680).

### REAGENTS

Reagents are stable until the date printed on the label assuming the container remains unopened at a storage temperature of 2-8 °C.

RUO LZ-SAA Calibrator EIKEN 1 mL x 6,

Ready to use.

### WARNINGS AND PRECAUTIONS

- For research use only.
- Do not use reagents that have passed their expiration date.
- Air bubbles on the surface of the product after being poured into sample cups can result in error measurement and should be removed.
- No human-derived ingredients are contained in this product, however, the risk of infection and the possible existence of other pathogens cannot be ruled out completely. Handle it as carefully as a patient's specimen. Treat it as a potential biohazard.
- Use the reagents as quickly as possible after they are opened. If they are to be stored, be sure to close the caps and store them using the prescribed method.
- Disposal of used reagents and containers should be handled as infectious waste in accordance with applicable regulations.
- \*7. The product contains Guanidinium chloride (CAS No.50-01-1) and ProClin 300 (CAS No.55965-84-9) which are classified by applicable 'The Classification, Labelling and Packaging (CLP) Regulation' as Skin and Eye Irritant and Acute toxicity by Oral. If the reagent contacts with eyes, mouth or skin, rinse it out with large volumes of running water, and perform other required first aid. If necessary, seek medical attention.



### Warning

H302	Harmful if swallowed.
H315	Causes skin irritation.
H317	May cause an allergic skin reaction.
H319	Causes serious eye irritation.
P280	Wear protective gloves / protective clothing / eye protection / face protection.
P333 +P313	If skin irritation or rash occurs: Get medical advice / attention.
P337 +P313	If eye irritation persists: Get medical advice / attention.
P362 +P364	Take off contaminated clothing and wash it before reuse.

### PROCEDURE

Create the calibration curve in the manner described by automatic analyzer manual.

### PRODUCT CODE, PRODUCT NAME & STORAGE

Product code	Product name	Contents	Storage
V-SZ12	RUO LZ-SAA Calibrator EIKEN	1 mL x 6	2-8 °C

### \*CONCENTRATION

Calibrator No.	Concentration (mg/L)
1	0.0
2	10.0
3	24.5
4	70.0
5	140.0
6	225.0



**EIKEN CHEMICAL CO., LTD.**

4-19-9 Taïto, Taïto-ku, Tokyo, 110-8408 JAPAN

\*Date of Revision: 2024-02-01

Date of Issue: 2021-07-01